


# IMPORTANT HANDLING INFORMATION

## For ibidi Heating System, Multi-Well Plates

# READ THIS BEFORE USE

Read this information sheet, in its entirety, **BEFORE** working with the ibidi Heating System for Multi-Well Plates.

The manual should also be read before using the instrument.

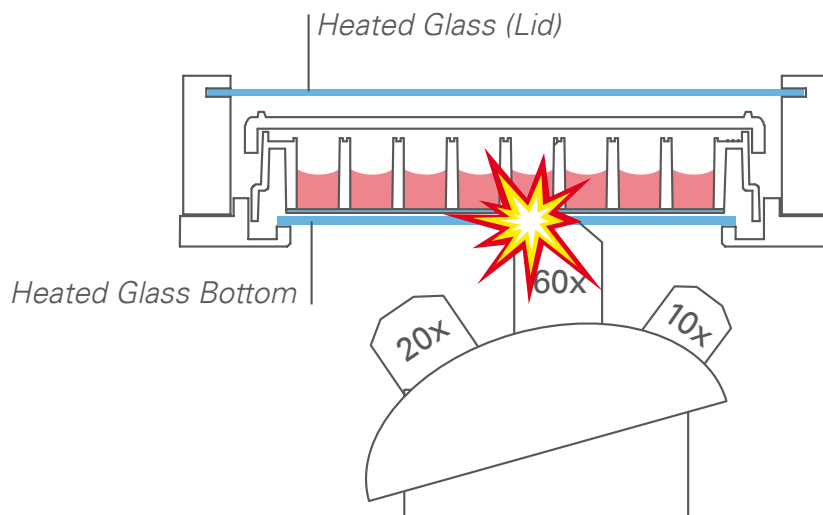

### CAUTION

The glass bottom of the heated plate may break due to a mechanical impact. Avoid contact between the glass bottom and the objective lens. Changing objective lenses may automatically lead to a collision between the objective lens and the glass bottom. Note: Never use an objective lens with a working distance smaller than 1 mm.

### CAUTION

If the glass plate of the heated lid or the plate itself breaks due to a mechanical impact, the shards can lead to injuries, if handled.